Prestained Protein Marker(10-250kDa)

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The 3-color Prestained Protein Marker Marker is a ready to use three-color protein standard with 11 pre-stained proteins covering a wide range of molecular weights from 10 to 250 kDa (10 to 230 kDa in Bis-Tris /MOPS buffer). Proteins are covalently coupled with a blue chromophore except for two reference bands (one **green** at 25 kDa and one **red** band at 70 kDa respectively) when separated on SDS-PAGE (Tris-Glycine buffer).

The Protein Marker is designed for monitoring protein separation during SDS-polyacrylamide gel electro-phoresis, verification of Western transfer efficiency on membranes, (PVDF, nylon, or nitrocellulose) and for approximating the size of proteins.

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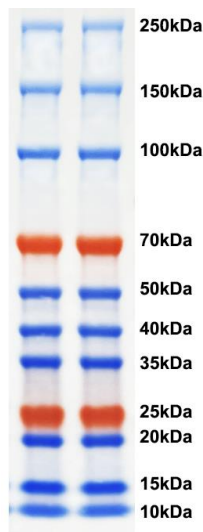
1. 产品已预混 loading buffer, **不需要加热、稀释或添加还原剂**。The marker is supplied in gel loading buffer and is ready to use. **Do NOT heat, dilute, or add reducing agents before loading.**
2. 加样前在室温放置数分钟, 解冻后, 轻轻摇匀, 以确保溶液混合均匀。Let it stand at room temperature for a few minutes before using, and after thawing, shake it gently to ensure that the solution is mixed evenly.

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